

## DSX Application Software Ver. 2.1.5

### Release Note

System Configuration	
OS	Microsoft Windows7 Professional 32bit/64bit SP1
	Japanese or English
CPU	Intel Xeon®E3-1230 Quad-Core3.2GHz or more
RAM	4GB or more
HDD	500G or more ※The capacity to save images, etc. by the DSX application software depends on the free space in HDD.
Graphic	Equivalent to NVIDIA Quadro 600 1GB or more.
Optical Drive	DVD drive (The offline software can be installed to CD drive.)
Monitor Resolution	1920 x 1080

**1. Changes from the previous version (Ver. 2.1.4)****1-1 Following phenomena were corrected.**

No	Function	TicketNo	Description
1	Particle analysis	#4929	Edit particles functions are disabled when strong/weak slider state changed during editing particles
2	User Calibration	#4992	Application crashes while exporting step measurement result to datasheet after changing user calibration X/Y values.
3	Wizard	#4984	Error that edges were not detected is displayed while executing switch image wizard and caliper wizard.
4		#4985	Application become unstable while editing caliper wizard node text box values using spin control.
5		#4989	Loading wizard data file of the OLS from disk is making application unstable.
6		#4991	Application become unstable while double click multiple times for drawing polygonal line for profile measurement.
7		#4923	Application hangs or is unstable when switch to reports happen from volume measurement where measurements are taken both in profile and histogram mode.
8		#4993	Application hangs when running CSV export wizard node and at same time respective CSV file is already open by operator.